



EPIX10KA

Based on e-mail exchange between Mikhail, Gabriel, Jack, Faisal, Philip.

For composite detectors see [EPIX10KA2M](#) and [EPIX10KAQUAD](#)

- [Pictures from Gabriel's presentation](#)
- [Configuration objects for epix10ka](#)
- [Gain coding](#)
 - [Data gain bits assignment](#)
 - [Pixel Mask](#)
- [Calibration rows](#)
- [Calibration files for dark](#)
- [Raw data](#)
- [Gain correction files](#)
- [Control bits table](#)
- [Gain correction formula](#)
- [Common mode correction](#)
- [Meeting with Gabriel](#)
- [Calibration Details](#)
- [Per panel calibration constants](#)
 - [File naming convention in panel calibration repository](#)
- [Constants merging for multi-panel detectors](#)
- [Pedestal Deployment Meeting Notes \(8/2/18\)](#)
- [List of known detectors on 2018-11-12](#)
- [Composite detectors](#)
- [References](#)

Pictures from Gabriel's presentation

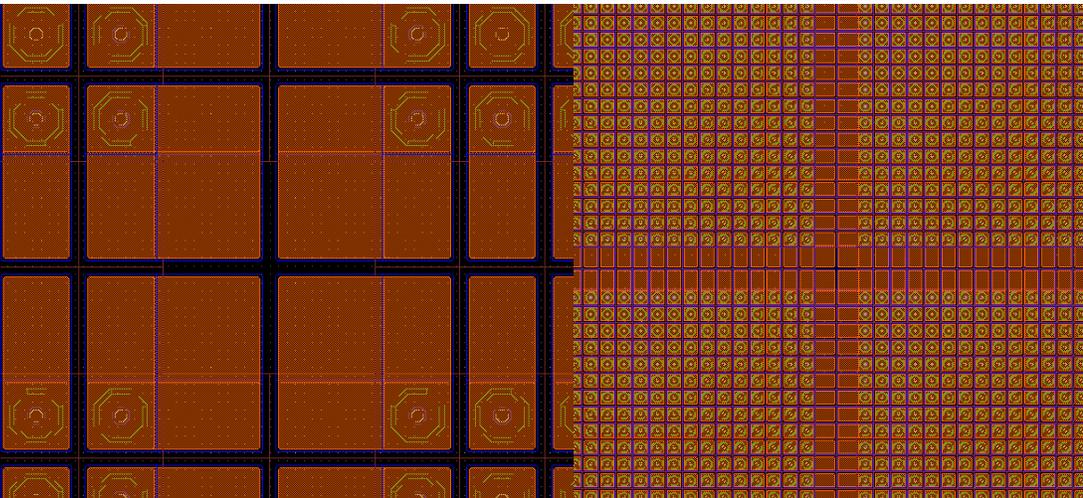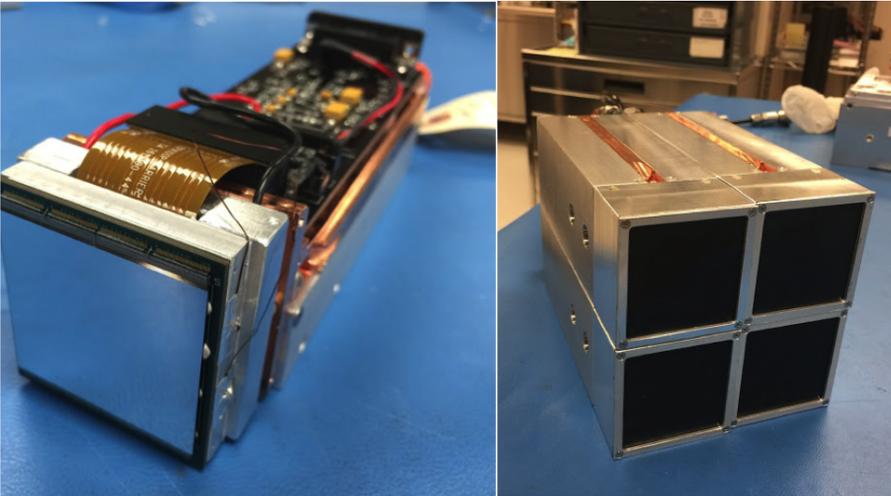

Chris Kenney:

- sensor regular pixel size 100x100 μm^2 ,
- shown on plot gap pixels 100x225 μm^2 and four central pixels 225x225 μm^2 .
- See plot for composite detectors in [EPIX10KA2M](#) and [EPIX10KAQUAD](#)

Configuration objects for epix10ka

Python psana has currently defined a few configuration objects for epix10ka:

psana.Epix.Config10kaV1		
co.acqToAsicR0Delay	co.asicMask	co.calibPixelConfigArray
co.numberOfCalibrationRows	co.scopeEnable	co.calibrationRowCountPerASIC
co.adcClkHalfT	co.asicPixelConfigArray	co.carrierId0
co.numberOfColumns	co.scopeTraceLength	co.carrierId1
co.adcPipelineDelay	co.asicPpbe	co.dacSetting
co.numberOfOfEnvironmentalRows	co.scopeTrigChan	co.digitalCardId0
co.adcPipelineDelay0	co.asicPpbeControl	co.digitalCardId1
co.numberOfPixelsPerAsicRow	co.scopeTrigEdge	co.enableAutomaticRunTrigger
co.adcPipelineDelay1	co.asicPpmat	co.environmentalRowCountPerASIC
co.numberOfReadableRows	co.scopeTrigHoldoff	co.epixRunTrigDelay
co.adcPipelineDelay2	co.asicPpmatControl	co.evrDaqCode
co.numberOfReadableRowsPerAsic	co.scopeTrigOffset	co.evrRunCode
co.adcPipelineDelay3	co.asicPPmatToReadout	co.evrRunTrigDelay
co.numberOfRows	co.SyncDelay	co.numberof125MhzTicksPerRunTrigger
co.adcReadsPerPixel	co.asicR0	co.numberOfAsics
co.numberOfRowsPerAsic	co.SyncMode	co.numberOfAsicsPerColumn
co.adcStreamMode	co.asicR0ClkControl	co.numberOfAsicsPerRow
co.prepulseR0Delay	co.SyncWidth	
co.analogCardId0	co.asicR0Control	
co.prepulseR0En	co.testPatternEnable	
co.analogCardId1	co.asicR0ToAsicAcq	
co.prepulseR0Width	co.TypeId	
co.asicAcq	co.asicR0Width	
co.R0Mode	co.usePgpEvr	
co.asicAcqControl	co.asicRoClk	
co.scopeADCsameplesToSkip	co.Version	
co.asicAcqLToPpmatL	co.asicRoClkHalfT	
co.scopeADCThreshold	co.version	
co.asicAcqWidth	co.asics	
co.scopeArmMode	co.asics_shape	
co.asicGR	co.baseClockFrequency	
co.scopeChanAwaveformSelect		
co.asicGRControl		
co.scopeChanBwaveformSelect		

psana.Epix.Asic10kaConfigV1 as co.asic 			
asic.atest	asic.FELmode	asic.RO_rst_en	asic.S2D_tcomp
asic.chipID	asic.Filter_DAC	asic.RowStart	asic.Sab_test
asic.ColumnStart	asic.Hrtest	asic.RowStop	asic.SLVDSbit
asic.ColumnStop	asic.is_en	asic.S2D	asic.tc
asic.CompEn_lowBit	asic.Monost	asic.S2D0_DAC	asic.test
asic.CompEn_topTwoBits	asic.Monost_Pulser	asic.S2D0_GR	asic.testBE
asic.CompEnOn	asic.OCB	asic.S2D0_tcDAC	asic.testLVDTransmitter
asic.CompTH_DAC	asic.Pbit	asic.S2D1_DAC	asic.TPS_DAC
asic.DelCCKreg	asic.PixelCB	asic.S2D1_GR	asic.TPS_GR
asic.DelEXEC	asic.pixelDummy	asic.S2D1_tcDAC	asic.TPS_MUX
asic.DM1	asic.PP_OCB_S2D	asic.S2D2_DAC	asic.TPS_tcDAC
asic.DM1en	asic.Preamp	asic.S2D2_GR	asic.TPS_tcomp
asic.DM2	asic.Pulser	asic.S2D2_tcDAC	asic.trbit
asic.DM2en	asic.Pulser_DAC	asic.S2D3_DAC	asic.Vld1_b
asic.emph_bc	asic.PulserR	asic.S2D3_GR	asic.VREF_DAC
asic.emph_bd	asic.PulserSync	asic.S2D3_tcDAC	asic.VrefLow
asic.fastPP_enable	asic.RO_Monost	asic.S2D_DAC_Bias	


```

#----|----#
# A2 | A1 #
#----|----#
# A3 | A0 #
#----|----#
# ||||| ||||| #
wirebonds

```

```

# r0 is row 176 in ASIC0 and ASIC3
# r1 is row 176 in ASIC1 and ASIC2

```

```

# r2 is row 177 (last) in ASIC0 and ASIC3
# r3 is row 177 (last) in ASIC1 and ASIC2

```

```

# when pixel (ga,g ,M,T) is xx0x then 176 is baseline and 177 is pixel max
# when pixel (ga,g ,M,T) is xx1x then 176 is pixel max and 177 is baseline

```

Calibration files for dark

2018-02-28 Philip:

"force switch" mode: the camera starts high then goes low - this simulates switching. This mode should be used for calibration of pedestals for switched mode.

I think by the way that we need to have two pedestal files in standard running:

H/M (dark)

H/M->L (forced)

In the case where we run the array in H/M/L without auto-ranging, we can just populate the relevant pedestals.

2018-02-28 Angelo:

The force switch mode should be ignored. It is a debugging mode that should not be used for calibration. Gabriel can explain how to do the calibration if needed

Raw data

epix10ka data available on 2018-07-05

```
Command: find_detector_runs MFX epix10ka
```

```


# mfx32516: nruns 3-377, MfxEndstation.0:Epix10ka.0, MfxEndstation.0:Epix10ka.1
# mfx36916: runs 1-84, MfxEndstation.0:Epix10ka.1

```

```

dataset      exp=mfx32516:run=377
calibDir:    /reg/d/psdm/MFX/mfx32516/calib
det.source   : Source("MfxEndstation.0:Epix10ka.0")
shape of ndarray: (352, 384)

```

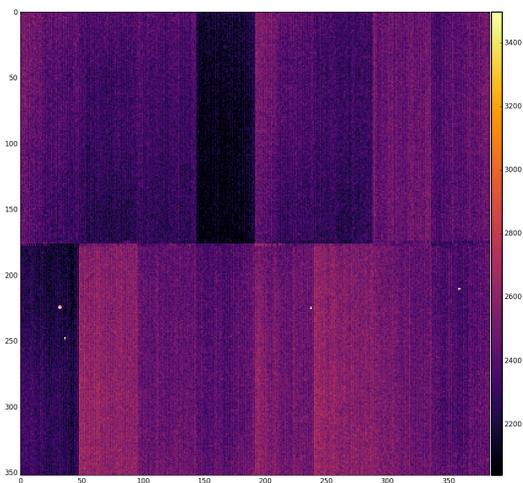

Gain correction files

On 2018-06-20 Gabriel generated gain and offset files in

gain correction files

```

Camera1-AHL-H-Gain.txt
Camera1-AHL-L-Gain.txt
Camera1-AML-L-Gain.txt
Camera1-AML-M-Gain.txt
Camera1-FH-H-Gain.txt
Camera1-FL-L-Gain.txt
Camera1-FM-M-Gain.txt
    
```

Archived in /reg/g/psdm/detector/gains/epix10k/2018-06-04-Camera1/

Gain and offset Mean and RMS

	Gain		Offset	
	Mean	RMS	Mean	RMS
FL-L	1.28	0.06	0	?
FM-M	42.43	2.00	2331	150
FH-H	128.49	6.17	2354	151
AML-L	1.26	0.06	2115	164
AML-M	42.52	1.92	2334	150
AHL-L	1.29	0.06	2119	190
AHL-H	128.72	6.03	2358	151

Control bits table

cbits - pixel control bit array

```

#-----
# data bit 14 is moved here 1/0 for H,M/L
# / trbit 1/0 for H/M
# V / bit3 1/0 for F/A
# V / bit2 1/0 for H,M/L
# V / M mask
# V / T test gain range index
# V / / in calib files
# V V
# x111xx =28 - FH_H 0
# x011xx =12 - FM_M 1
# xx10xx = 8 - FL_L 2
# 0100xx =16 - AHL_H 3
# 0000xx = 0 - AML_M 4
# 1100xx =48 - AHL_L 5
# 1000xx =32 - AML_L 6
# 1101xx =49 - AHL_FL 7
# 1001xx =33 - AML_FL 8
#-----
    
```

Index	Gain mode	Data	config	asicPixelConfigArray			
		bit14	trbit	bit3: ga	bit2: g	bit1: Mask	bit0: Test
0	FH_H	x	1	1	1		
1	FM_M	x	0	1	1		
2	FL_L	x	x	1	0		
3	AHL_H	0	1	0	0		

4	AML_M	0	0	0	0		
5	AHL_L	1	1	0	0		
6	AML_L	1	0	0	0		
7	AHL Forced L	1	1	0	1		
8	AML Forced L	1	0	0	1		
pixel control bit array		$1 \ll 5 = 32$	$1 \ll 4 = 16$	$1 \ll 3 = 8$	$1 \ll 2 = 4$	Not used	Not used

Gain correction formula

corrected = (raw - pedestals)/gain

Common mode correction

Added on 2020-06-04.

The value (raw - pedestals) can be corrected for **common mode** effect.

First image below shows one epix10ka segment non-corrected image of shape (352, 384). Vertical stripes presumably arising from this hardware effect.

The same median algorithm is applied as for Jungfrau, but separately for up- and down- part of the segment with parameters

cmpars=(**<algorithm-id>**, **<mode>**, **<maximal-correction>**), where

- algorithm-id is not used
- mode=0 - no correction, bit1=1 - in rows (is not useful for epix10ka), bit2=1 - in columns (the best for epix10ka)
- maximal-correction - maximal allowed correction in ADU

NOTE: currently mask on gain mode is not applied - all pixels are used to evaluate and correct for offset.


Examples of the command to get corrected image:

```
nda_cdata = det.calib(evt, cmpars=None) # default - cm pars from calib directory/system
```

```
nda_cdata = det.calib(evt, cmpars=(0, 0, 100)) # use cmpars, do not apply cm correction
```

```
nda_cdata = det.calib(evt, cmpars=(0, 2, 100)) # use cmpars, apply cm correction in columns
```

Images before and after common mode correction:

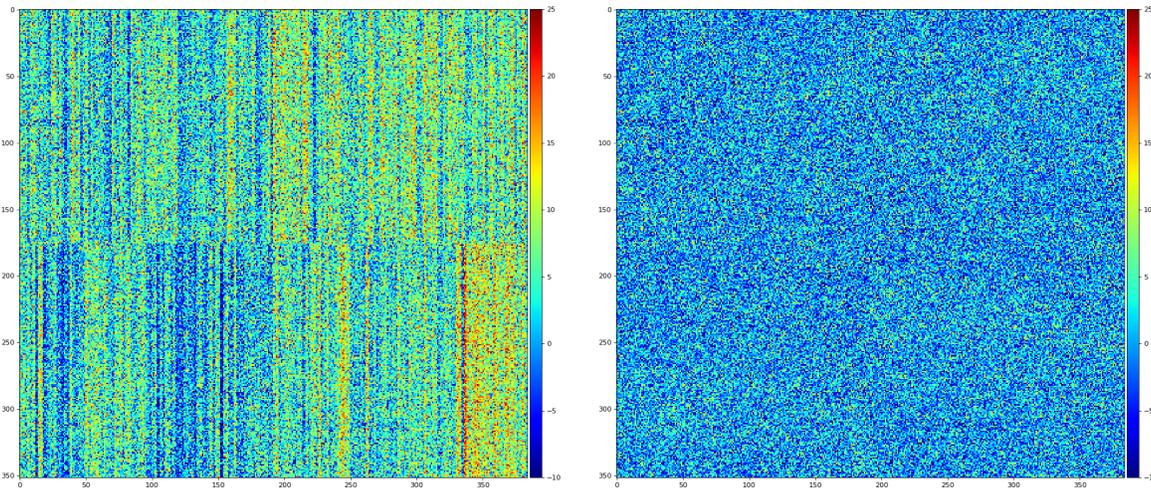

Meeting with Gabriel

On 2018-07-06 15:13 meeting with Gabriel.

Summary:

- control bit table is correct
- formula should include pedestals in stead of offset

- pedestals need to be generated and saved as txt file of shape (7, 1, 352, 384) under calib directory, for example: /reg/d/psdm/MFX/mfxx32516/calib/Epix10ka::CalibV1/MfxEndstation.0/Epix10ka.0/pedestals/0-end.data
- Gabriel on pedestal calibration:
 - offsets and gain are calibrated not very often (say, once per month...)
 - pedestals should be calibrated pretty often (once per hour...)
 - pedestal calibration is implemented in AMI, using offsets,
 - script works on dark runs accumulated for 5 gain modes
 - + 2 gain modes are evaluated using combination of other 2 and offsets
 - pedestal calibration procedure needs to be completed.

Calibration Details

See this page: <https://confluence.slac.stanford.edu/display/ppareg/ePix10K+Calibration>

Per panel calibration constants

Algorithm for xtc data with charge injection processing is developed by Gabriel and is wrapped in psana environment as CLI

Detector/app/epix10ka_offset_calibration

Similarly, dark run processing script is

Detector/app/epix10ka_pedestals_calibration

These commands will be available in releases > ana-1.3.58. See description using option "-h", e.g.

```
command> epix10ka_offset_calibration -h
```

```
command> epix10ka_pedestals_calibration -h
```

File naming convention in panel calibration repository

/reg/g/psdm/detector/gains/epix10k/panels/ - default repository <dir-repo>, can be re-directed for test by option "-o <dir-name>".

<dir-repo>/3791847426-0170080513-1879048214-0191557724-0003673288-2996154369-0218103833/ - directory for particular panel includes its long id,

Sub-directories for each panel:

- offset
- pedestals
- plots
- work
- gain # content will be provided by Philip

<dir-repo>/aliases.txt - file with aliases for long panel ids,

<dir-repo>/panels/logs/ - directory for log files,

Panel calibration file name: epix10ka_0001_20180514120622_mfxx32516_r1021_pedestals_AHL-H.dat, where

- epix10ka - detector type
- 0001 - panel id alias as specified in /reg/g/psdm/detector/gains/epix10k/panels/aliases.txt
- 20180514120622 - run start time stamp in format '%Y%m%d%H%M%S'
- mfxx32516_r1021 - experiment and run for production of this constants
- pedestals - constants type
- AHL-H - gain range (mode)

File with already processed arrays is saved as

```
<dir-repo>/<panel-id>/work/epix10ka_0001_20180514120622_mfxx32516_r1021_sp02-df.npz
```

with extra fields in name sp02 for charge injection spacing number of pixels.

Constants merging for multi-panel detectors

```
command> epix10ka_deploy_constants -h
```

shows description of command line options. This command with sufficient number of parameters (at least experiment, run, and detector name)

- combines gain mode and panel calibration files in arrays of shape (7, <number-of-panels>, 352, 384)
- deploys calibration files in the calib directory, e.g.

```
/reg/d/psdm/MFX/mfxx32516/calib/Epix10ka::CalibV1/MfxEndstation.0:Epix10ka.0/pixel_gain/0-end.data
```

```
/reg/d/psdm/MFX/mfxx32516/calib/Epix10ka::CalibV1/MfxEndstation.0:Epix10ka.0/pedestals/0-end.data
```

Pedestal Deployment Meeting Notes (8/2/18)

- script that is use to drive the daq with hutch python: `/reg/g/pcds/pyps/mfx/dev/mfx/experiments/mfxls3416.py`
- for epix, 5 gain ranges (3 fixed, 2 auto) generate 7 pedestals
- maybe have "all range" option, and fast "selected ranges" option.
- maybe "all range" at beginning of expt, but "selected ranges" during expt.
- jason thinks medium/low is the default
- in the "selected ranges" mode makepedes will reuse the current numbers for ranges that weren't selected.
- multi-panel epix10k should be the same as jungfrau (might need to massage the daq appropriately)
- mikhail/gabriel: data processing call in makepedes script
- silke/jason/dan: hutch python, makepedes wrapper
- `/reg/g/pcds/engineering_tools/R1.2.9/scripts/makepedes` calls `makepedes_psana` calls mikhail's stuff (`jungfrau_ndarr_dark_proc`)

List of known detectors on 2018-11-12

```
command> find_detector_names
```

```
#== type 0022 epix10ka
      'DetLab.0:Epix10ka.0': 'epix10ka_0000',\
      'MecTargetChamber.0:Epix10ka.0': 'epix10ka_0001',\
      'MecTargetChamber.0:Epix10ka.1': 'epix10ka_0002',\
      'MfxEndstation.0:Epix10ka.0': 'epix10ka_0003',\
      'MfxEndstation.0:Epix10ka.1': 'epix10ka_0004',\
      'MfxEndstation.0:Epix10ka.2': 'epix10ka_0005',\
```

Composite detectors

[EPIX10KA2M](#) and [EPIX10KAQUAD](#) - composite detectors made of epix10ka panels.

References

- <https://github.com/slaclab/cac/tree/master/psana/epix10ka> - scripts from Faisal
- https://pswww.slac.stanford.edu/apps/portal/index.php?exper_id=1107 - under Workflow -> Batch Processing - batch processing
- `/reg/d/psdm/MFX/mfxx32516/results/abunimeh/2` - experimental and jupyter notebook scripts
- [ePix10K Calibration](#), by Gabriel Blaj
- [EPIX10KA2M and EPIX10KAQUAD](#)
- [Jungfrau and Epix10ka Calibration](#)
- 2021-09-28 learn about existence of presentation [2019-10-02-ePix10K-500um-Sensor-Geometry.pdf](#)